

Testdigest

Name:

Date:

Continue from Experiment
(Date)

(Name)

Project Name:

For one reaction you need: For Mastermix: Number of
samples+2extra

4µl H₂O

1µl Buffer, NEB4

1µl BSA (10x)

0,5 µEnzym 1

0,5 µEnzym 2

3 µl DNA

10 µl total volume

Give 3 µl of DNA in an eppi and add 7µl of the mastermix.
Incubate for about 1h at 37°C.

Add 1 µl Loading dye buffer and load the gel.

Take a picture of the gel, print picture and label the lanes!